


From hot to cold

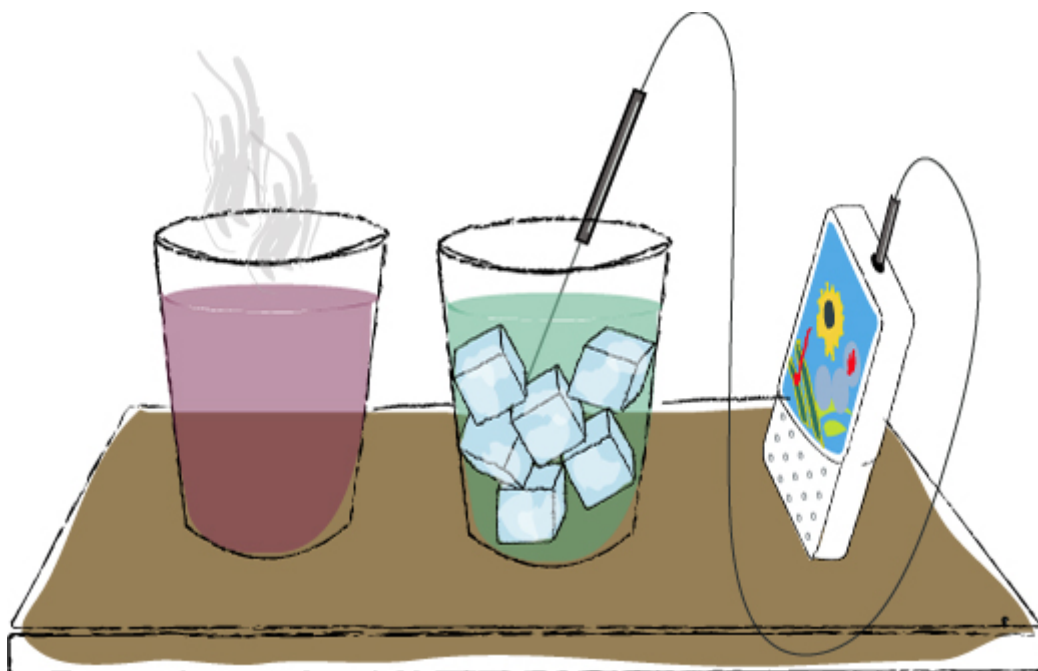

We're going to...


Explore the use of a temperature sensor and the way it responds to changes in temperature

Getting started

- ✚ Connect the temperature sensor to IN - A on EcoLog XL
- ✚ Click **Display Workbook**  on the main toolbar.
- ✚ Click **Lets Go...** then click **From hot to cold**
- ✚ To start the experiment click **Launch**  on the lower toolbar

To show or hide the workbook click **Workbook**  on the main toolbar


Let's go


1. Click **Run**  on the upper toolbar to begin
2. Place the temperature sensor in the hot water
3. Watch the changes on each display
4. Place the temperature sensor in the cold water
5. Again, watch the changes on each display



What did we learn?

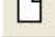

Click **New**  on the lower toolbar to open a new text window and type in your answers to the following questions

(don't forget to save your answers - click **Save**  on the lower toolbar):

1. Which display best shows the changes?
2. How hot is the hot water?
3. How cold is the cold water?

